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CONTINUATION

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Description in the manual is based on the newest products. Owing to improvement or other changes, contents of manual may differ from practical situation. Our company will reserve the right of recension at any moment. Please kindly forgive not to notify the revised place one by one.

The copyright of this manual belongs to Wenzhou Sundoo Instruments Co., Ltd. Any companies or individuals have no right to copy or plagiarize part or whole contents of the manual without permission of our company.

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MANUFACTURER: WENZHOU SUNDOO INSTRUMENTS CO.,LTD

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Thank you for your purchasing SVM-212 Stereo Microscope.

SVM-212 Stereo Microscope is researched and designed for quality control of modern manufacture. It can provide high clear image by adopting unique technology of eliminating interferential beam and wide range field, and Hiquality SONY CCD camera. Moreover the item owns up and down double illumination, and can be used in various environment. The video of microscope can be directly output to view, which helps inspector to avoid optic tiredness and reading from scale. It also realizes measurement of length,radius, angle, etc. In some aspect, the microscope can replace profile projector, industrial microscope and image survey meter with easy operation. This product is widely used in fields of electron, spring, hardware, mould, powder metallurgy, precision instrument, jewellery, clock, fiber, quality control, biology, medicine, biochemical study, and so on.

Packing List

Main body		Quantity
Upper and background illuminations		1
Lens base		1
Flat-field achromatic object lens	16×	1
Flat-fielt ocular	20×	2
Carry object platform		1
CCD Photographic system		1
User's manual		1
Package list		1
Certificate		1

(2) Display + Computer + Video Card + Measuring Software

It can realize micro measuring function by measuring software, which can measure width, radii/diameter, distance between two lines, angle etc. Connecting Method:

Please insert Video Card into PCI socket of computer, and install Video Card's drive program and measuring program. Then insert one head of AV cable into CCD AV output connector, and connect another head into input connector of Video Card.

(3) Moving Platform

Adding moving platform on work table will make measuring objects moving more smoothly. It attaches a scale with 0.02mm resolution on moving platform.

Maintains

- 1. When the microscope is not used, please screw ocular protector well, and cover dustproof to prevent dirty, then put in dry place. We advise to stock binocular (trinocular) base and ocular into dampproof container. and place drier to prevent mould.
- 2. When the microscope is cleaned, don't touch the optical accessories by hands. Please use soft cotton or brush dipped with pure alcohol to clean dirt, obstinate dirt and fingerprint, grease etc.

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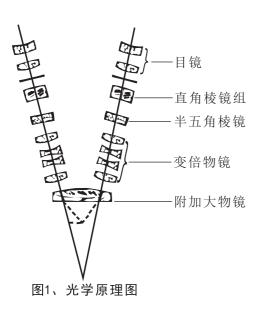
lower magnification can provide wider field of view and depth of scene, which is suitable to find object. We advise to find inspected object by lower magnification at first, then turn Magnification Adjustable Handwheel to get higher magnification for inspecting the details of object.

(4) When magnification is changed from lower to higher, please turning Adjustable Handwheel slightly if the image is little unclear. If image is not clear during changing magnification, please check whether the optic angle is adjusted correctly.

6. Application of Lens Cone

When photography or colorful television is equipped on lens cone, please pull out pull rod of microscope base to make image on the straight cylinder. Loosen fixture screw on straight cylinder, and turn adjustable ring to change length. Then the image in straight cylinder will be same to slant cylinder (it is adjusted well when out of factory).

7. Changing Light Bulb

Please put off the power before changing light bulb. After the lamp is cold, remove Slant Illumination Cover, take off halogen bulb and change it.

Opening window on the bottom base can change halogen or fluoresce bulb.

8. Application of Accessorial Ocular

The accessorial ocular is optional to match with microscope if requiring longer working distance or larger magnification. It will have longer working distance by 0.5x and 0.75x, and will own larger magnification by 1.5x and 2x. When using, just screw accessorial ocular into binocular (trinocular) base.

Optional Attachment

(1) AV/VGA, XGA-transfer card + CRT display

The XGA-transfer card can provide hi-definition image of 1024x768x75 Hz, display on CRT without flicker but with bright and flamboyant image. Connecting Step:

a. Attach the power, connect the CRT's signal connector with the transfer card's connector.

b.Inert matched MMI cable into connector on transfer card, then insert one terminal of AV cable into output connector of CCD camera, and insert another terminal into yellow connector of MMI cable.

SUNDOO® 世 仪器

Operation

1. Installation

Please take binocular (trinocular) ocular base and main body out of packing box. Lock Fixture Handwheel tightly to prevent Up and Down Pedestal falling down. Please fix the binocular (trinocular) base on main body by fixture screw, and take Lens Protector off.

2. Check Input Voltage

Please check whether voltage of power supply is according with rate input voltage of microscope, which is symbolized on the mark of back or side of the microscope base. If they are identical, put through the power, and the socket must have the earth wire.

3. Select Illumination

Please select slant illumination or transparent illumination according to inspecting requirements. The two illumination mode can be applied simultaneously or separately. The brightness of slant lamp and transparent lamp can be adjusted by Brightness Adjustable Handwheel. If transparent lamp is fluorescent light, the brightness can not be adjusted. The angel of slant illumination is adjustable to ensure enough bright of theinspected sample.

4. Adjust Pupil-gap

Turn Prism Cover to adjust horizontal distance between two oculars according to inspector's pupil gap. When eyes overlap beam through oculars, inspector can see left and right image overlapping.

5. Adjusting Focus

- (1) When inspector's sight is normal. Turn Optic Angle Adjustable Ring to zero. If the inspector is hyperopia or short sight, turn Optic Angel Adjustable Ring to corresponding place. For example, if left eye sight is myopia 200 degree, the optic ring should be turned to scale of . If the slight is hyperopia 200 degree, please turn the optic ring to scale of . Up and Down Pedestal Adjustable Handwheel
- (2) Adjust focus by turning Adjustable Handwheel, or changing height of Up and Down Pedestal.
- (3) Turning Magnification Adjustable Handwheel can provide continuous adjustable magnification. The magnification of object lens is 0.7x-4.5x. The

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2. Check Voltage

- (1)The rate input voltage is signed on the pedestal of microscope. The local power supply must be according with the rate input voltage. Otherwise, please ask the experience specialist to open the cover on the pedestal, turn the switch which has symbol of 110V, 220V, and select the voltage same as the power supply, then fix the cover on the pedestal. It will damage the microscope if it is not operated with rate voltage.
- (2) The power socket should have earth cable.

3. Prevent Fire or Scald

- (1) The bulb on microscope is in high temperature during lighting. Pay more attention to not touch the bulb to prevent scald.
- (2) When changing light bulb, it should be proceeding after bulb is completely cold.
- (3)Do not take tinder closing to bulb to avoid fire.

4. Change Light Bulb

- (1)Please cut off power before changing bulb to prevent electric shock or damage of equipment.
- (2) The microscope only adopts halogen bulb or fluorescent lamp.
- (3)Please wear gloves or pack the bulb with cover when changing light bulb, and do not touch the glass of the bulb by hand. After finished, please sweep finger mark and dirty on the bulb with clean cloth dipping in pure alcohol. Otherwise, the bulb will be easy to crack or brightness will be reduced.

5. Carry and Stock

- (1) The power must be cut off before moving the microscope.
- (2) Operation Environment:

Temperature: $0^{\circ}C-40^{\circ}C$.

Max Relative Humidity: 85%.

- (3)Don't put microscope under sunshine or lamp. bright environment will effect image quailiy.
- (4)Don't put microscope in dust environment, and it should be covered by dustproof when not use.
- (5)Please put microscope in place without vibration.

Specification

Model	SVM-212 Digital Stereo Microscope		
Image Device	Optic System1	1/3"CCD Sony	
Magnification	15×~90×Changing Magnification Continuously	50×~320×Changing Magnification Continuously	
Accuracy for Max Magnification	2 μ	2 μ	
Definition for Max Magnification	0.5 μ	0.5 μ	
Max Sight Field(mm)	20×20	6.5×4.9	
Display	Eyes Inspect	Monitor, Computer, Television, LCD	
Illumination	Upper & Background Illumination System		

Adjustable Range of Pupil-gap: 55~75mm; Adjusting focus range of handwheel: 60mm; Adjustable range of fluctuation: 110mm; Adjustable range of Optic Angel: ±5refraction; Illumination: Input voltage: 110V or 220V;

Output: Slant illumination: 12V 10W halogen bulb;

Transparent illumination: 12V 10W halogen bulb (Provide 5W

halogen bulb according to requirements)

Cautions

1. Don't Disassemble Microscope

To reduce the risk of electric shock and equipment damage, please do not remove accessories, which are not described to be removed in the manual. If malfunction, please refer service of professional maintenance man.

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
Picture 1 shows the optical theory of stereo microscope. It is composed with two completely identical beam paths, which are in an angel. The upper illumination and transparent illumination are offered with the microscope. The magnified and positive stereo image is inspected through the microscope. By optional purchasing Accessional Object Lens, it can obtain higher magnification or working distance. Most importantly, it has added one viewfinder and right-angle prism to make image on straight cyli

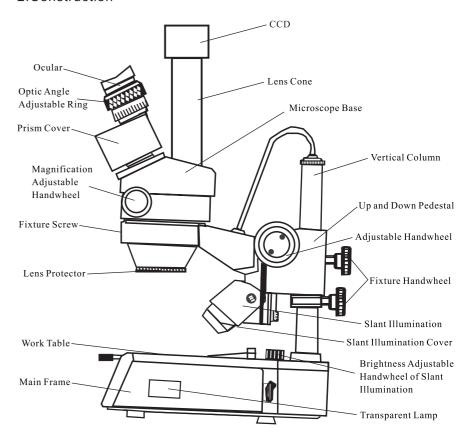2. Construction

Picture 2 Construction of SVM-212

Summary

1. Function

SVM-212 Digital Stereo Microscope is binocular and trinocular microscope changing magnification continuously. It can zoom the small object, and form stereo image. The microscope is widely used in electric industry, inspection, repair of small precise accessories, jewellery check and biological research. It can be applied with photography equipment or colorful micro-television.

2. Character

- (1)SONY CCD with high definition
- (2)Large work table
- (3) High work distance
- (4) Double lamps (background lamp, bottom lamp)
- (5) Video output (Realize measuring function by matching measuring software)

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